


Application Note

Phosphine & Arsine

Baseline - MOCON, Inc.™

1. Phosphine
2. Arsine

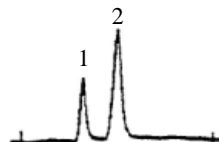

Chromatogram

Analysis

Detector: PID

Analysis Time: 2 min


Column: Proprietary

Oven Temperature: 65°C

Carrier: Nitrogen, 45cc/min typical

Represented by:

*Baseline, the reference point
from which all things are measured.*

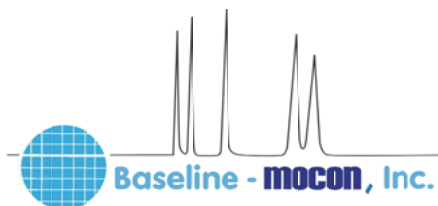

19661 Highway 36, P.O. Box 649, Lyons, CO 80540
In the continental United States, phone 800.321.4665, or fax 800.848.6464, toll free.
Worldwide, phone 303.823.6661 or fax 303.823.5151
• URL: www.baselineindustries.com • E-mail: sales@baselineindustries.com